



## PARAFFIN TRIMMER BHM1Y1 (BHTP-605)

# PARAFFIN TRIMMER BHM1Y1

## PARAFFIN TRIMMER



The grooved heating surface channels melted wax to a removable collection tray. The collection tray connects magnetically to the base of the main unit. Disposable plastic liners are included to insert into the collection tray for easy disposal of runoff wax. This unit now includes a digital control panel for personalized settings.

1. Remove the wax trimmer from its packaging and place it on a stable workstation. We recommend that you retain the carton and packaging.
2. The wax trimmer comes with a detachable AC power cord. Connect the cord to the port located on the back of the unit. When ready to use, plug the cord into a power source.
3. Attach the magnetic paraffin collection tray at the bottom of the unit and insert a plastic

## SPECIFICATIONS

|                                 |   |
|---------------------------------|---|
| Model                           | BHM1Y1  |
| Old Model                       | BHTP-605  |
| Power Switch & Display          | Flip the Power Switch to ON, digital display shows current temperature                |
| Temperature Setting             | SET Key to switch to temperature mode, adjust with side buttons, confirm with SET Key |
| Suggested Operating Temperature | 80-90°C (Do not exceed 100°C)   |
| Temperature Control Range       | Ambient temperature ~ 110°C   |
| Heating Indicator               | Red LED Light indicates heating   |
| Heating Time                    | Up to 10 minutes to reach desired temperature   |
| Safety Instructions             | Do not allow non-cassette items to touch surface, keep sleeves/clothing/jewelry away  |
| Cassette Handling               | Cassettes must be flat, pull downward to remove excess paraffin                       |
| Net Weight                      | 3 kg  |
| Instrument Dimensions           | 225*210*140 mm  |
| Package Info                    |   |
| --Packing                       | Wooden Case   |
| --Gross Weight                  | 4 kg  |
| --Package Dimensions            | 370*210*260 mm  |
| Alt Name                        | Paraffin Trimmer  |

## FEATURES

The grooved heating surface channels melted wax to a removable collection tray. The collection tray connects magnetically to the base of the main unit. Disposable plastic liners are included to insert into the collection tray for easy disposal of runoff wax. This unit now includes a digital control panel for personalized settings.

1. Remove the wax trimmer from its packaging and place it on a stable workstation. We recommend that you retain the carton and packaging.
2. The wax trimmer comes with a detachable AC power cord. Connect the cord to the port located on the back of the unit. When ready to use, plug the cord into a power source.
3. Attach the magnetic paraffin collection tray at the bottom of the unit and insert a plastic collection tray insert. The brim of the plastic liner can be folded down along the side of the tray that sits against the unit. The magnet will hold the collection tray in place.
4. Flip the power switch to the ON position to begin use and adjust temperature. (See Operation section for instructions).



**Biolab Scientific Ltd.**

Trillium Executive Center, East Tower, 675 Cochrane Dr, Markham, Ontario L3R 0B8, Canada  
Email: [info@biolabscientific.com](mailto:info@biolabscientific.com) | Website: [www.biolabscientific.com](http://www.biolabscientific.com)