



## VORTEX MIXER

BFA1BL1

# INDEX

1. Introduction	2
2. Product Features	2
3. Operation Instruction	3
4. Maintenance	4
5. Safety Warning	4
6. Parameters	5
7. Packing List	5

# 1. Introduction

The Touch Fairy Vortexer focus on touch-to-mix (short mix) for a variety of tubes / centrifugal tubes. Its compact design makes it very attractive and welcome, while its touch-to-mix operation makes it very easy handling and quite convenient. Touch fairy vortexer becomes an essential experimental instrument with its prominent mixing effect.

## 2. Product Features

1. The maximum diameter of the mixed tube is up to 30mm. Suit for mixing tubes / centrifugal tubes with prominent mixing effect.
2. Outer shell and touch concave are made of plastic which is resistance to chemical attack. The thick base makes operation very steady and reliable.
3. 12V low voltage power adapter built in, very smart and portable.

### Performance Test Statement

Name	Touch Fairy Vortexer	Type	BFA1BL1
Test Date		Production No.	MU-E48-0040
No.	Test Content	Standard	Test Results
1	Speed	4000±100 rpm	<input type="checkbox"/> Qualified
2	Basic Function	Valid	<input type="checkbox"/> Qualified
3	Appearance Demand	Valid	<input type="checkbox"/> Qualified
4	Appearance Sign	Valid	<input type="checkbox"/> Qualified
5	Continuous Work Test	72 hours no irregular	<input type="checkbox"/> Qualified
Test Result			
Tester:		Confirmer:	

Table 1

### 3. Operation Instruction

Put the instrument on a steady and horizontal bench top, connect the Power. Hold the tube to press the concave rubber , the instrument begins operating.

The vortex speed of BFA1BL1 is fixed.

### 4. Maintenance

Use mull with non-corrosive cleansing fluid (PH=7±1) to clean surface of the instrument. Operate the instrument after it dry out. Do not dip the instrument into fluid or water it.

### 5. Safety Warning

1. Do not use combustible, explosive and corrosive goods to the instrument.
2. The instrument should be put on a stable and firm working stable.

### 6. Parameters

Dimension: 110x110x86 mm	Net Weight: 1.1Kg
Power Input: AC90-230V, 50/60Hz	Speed: 4000rpm
Power: 12W	

Table 2

### 7. Packing List

	Name	Type	Quantity	Remarks
1	Touch Fairy Vortexer	BFA1BL1	1	
2	Adapter	12V/1A	1	
3	Performance Test Statement		1	
4	Operation Manual		1	
5	Warranty Card		1	
Charger: (Sign/Stamp)		Packing Date:		



**Biolab Scientific Ltd.**

Trillium Executive Center, East Tower, 675 Cochrane Dr, Markham, Ontario L3R 0B8, Canada  
Email: [info@biolabscientific.com](mailto:info@biolabscientific.com) | Website: [www.biolabscientific.com](http://www.biolabscientific.com)