



BIOLOGICAL MICROSCOPE

BMIC-205-A

INDEX

1. User Notice	2
1.1. Safety Note	2
1.2. Maintenance and Care	3
2. Components Name	4
3. Assembly	5
3.1 Assembly Diagram	5
3.2 Assembly procedure	6
4. Adjustment and operation	10
4.1 Adjustment set diagram	11
4.2 Adjustment and Operation	12
5. Technical Specifications	17
5.1 Main Specifications	17
5.2 Eyepiece and objectives	18
6. Outfit	19
7. Troubleshooting Guide	21
7.1 Optical System	21
7.2 Mechanical System	22
7.3 Electrical System	23

1. User Notice

1.1. Safety Note

1. Open the box carefully to avoid the accessories, like lenses, dropping to the ground or being damaged.
2. Do keep the instrument out of direct sunlight, high temperature or humidity, dusty and easy shaking environment. Make sure the stage is flat, horizontal and firm enough.
3. When moving the microscope, carefully carry it with the handle and the base.
4. Make sure the instrument is earthed, to avoid lighting strikes.
5. For safety, be sure the main switch is in "O" (off) state and cut off the power supply before replacing the bulb or the fuse. If you replace the bulb during use or right after use, allow the lamp bulb and the lamp house to cool completely before touching.
(Designated bulb:1W S-LED or 6V20W Halogen Lamp)
6. Check the input voltage: be sure the input voltage which is signed in the back of the microscope is consistent with the power supply voltage, or it will bring a serious damage to the instrument.
7. Always use the power cord provided by Novel.
8. The electrical equipment of the microscope should be discard as electronic waste.

1.2. Maintenance and Care

1. All the lenses have been adjusted properly; do not dismount them by yourself please.
2. The nosepiece and coarse and fine focusing parts are so delicate that it is forbidden to disassemble them carelessly by yourself.
3. Keep the instrument clean, and do not pollute the optical element when wiping away the dust on the instrument.
4. The contaminations on the prism, like fingerprints and oil smudges, could be gently wiped with a piece of soft cloth or tissue paper, gauze which has been immersed in pure alcohol or ether. (Note that the alcohol and ether are highly flammable, do keep them away from the fire or potential sources of electrical sparks, and use them in a drafty room as possible as you can.)
5. Do not attempt to use organic solvents to clean the microscope components other than the glass components. To clean them, use a lint-free, soft cloth slightly moistened with a diluted neutral detergent.
6. When using, if the microscope is splashed by liquid, cut off the power at once, and wipe away the splash.
7. Do not disassemble any parts of the microscope, as this will affect the function or reduce the performance of the microscope.
8. Place the instrument in a cool, dry position. When not using the microscope, keep it covered with a dust cover. Make sure the lamp socket is cool before covering the microscope.

2. Components Name

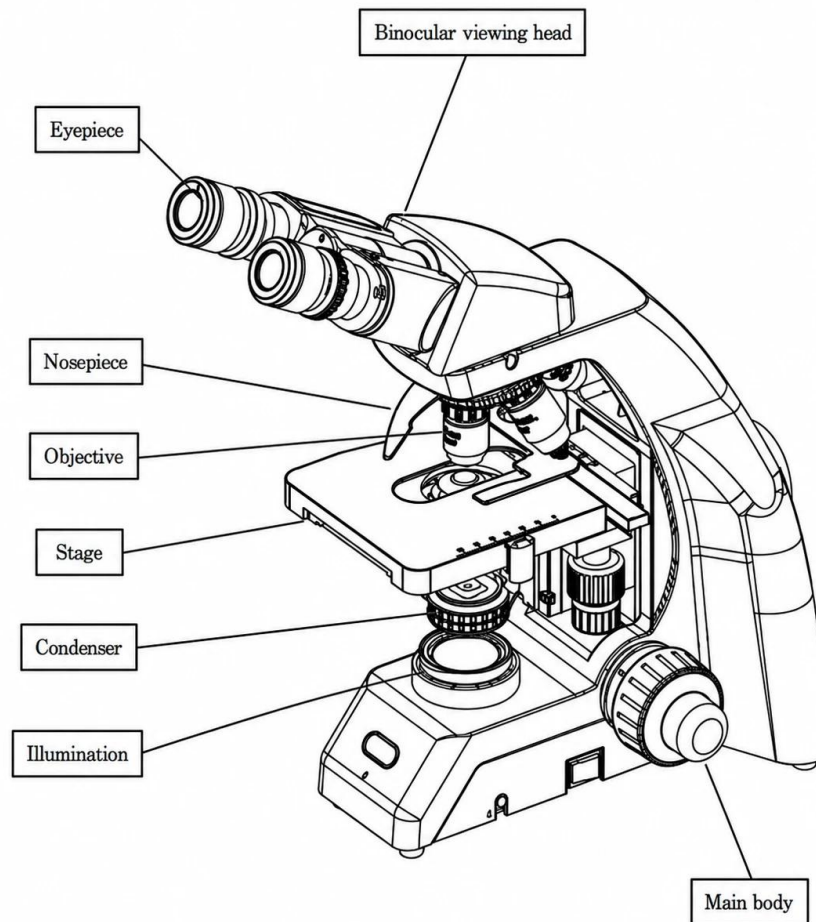


Figure 1

3. Assembly

3.1 Assembly Diagram

The following figure shows the installation sequence of the components. The number in the figure shows the assembly steps.

Before installing, be sure every components is clean, do not score any parts or glass surface.

Keep well with the hexagon wrench provided. When changing the components, you will need it again.

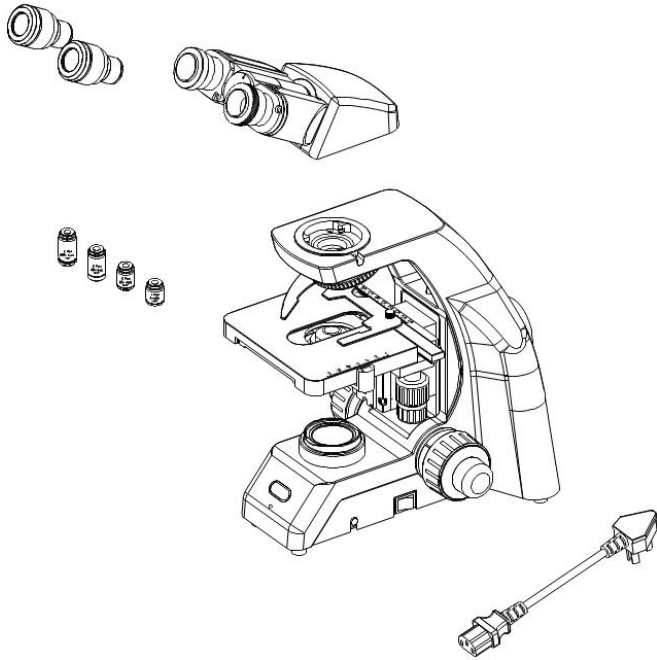


Figure 2

3.2 Assembly procedure

3.2.1 Installing binocular viewing head

Insert the binocular viewing head into the microscope head and turn it to a proper position, then tight it.

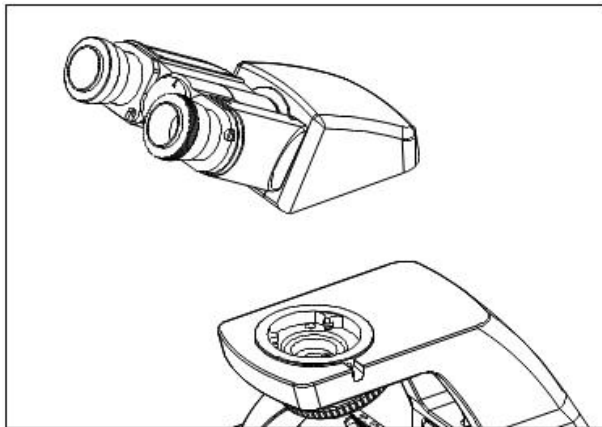


Figure 3

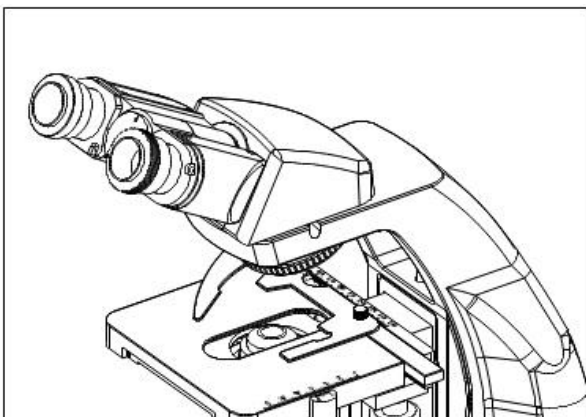


Figure 4

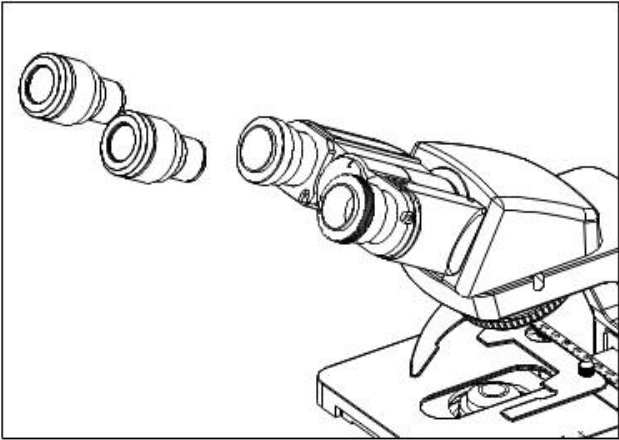


Figure 5

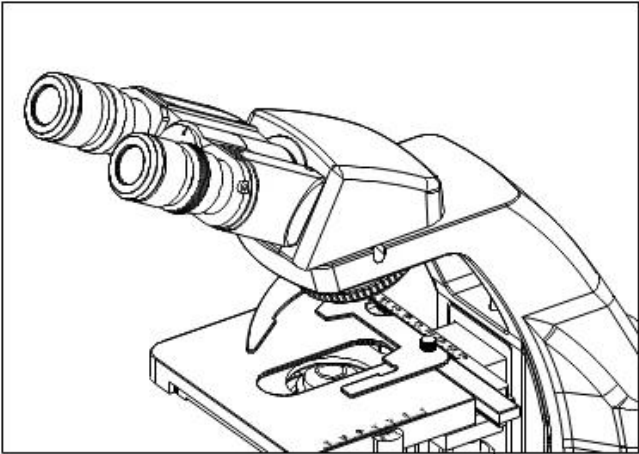


Figure 6

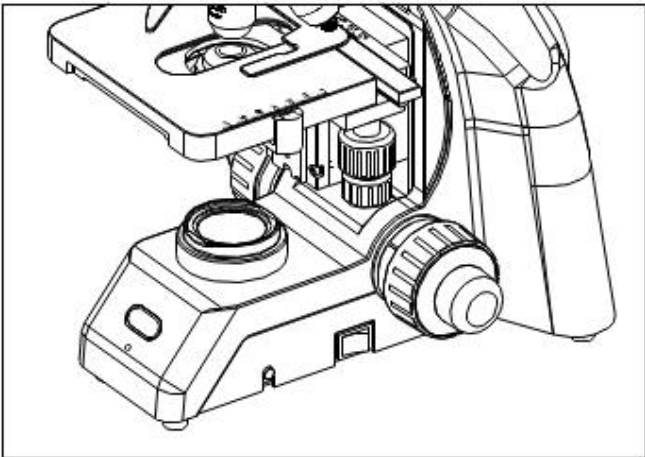


Figure 7

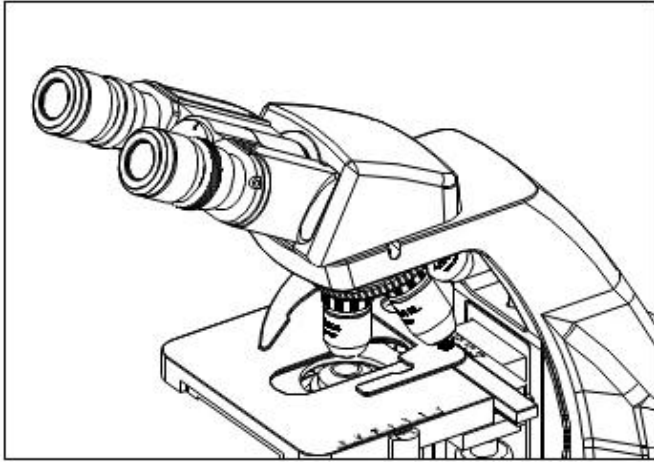


Figure 7

3.2.3 Installing the objective

1. Adjusting the coarse focus knob until the support device of the mechanical stage reaches its low limit position.
2. Screw the lowest magnification objective into the nosepiece from the left or the right side, then revolve the nosepiece clockwise and mount other objectives by the sequence of low to high magnification

Installing objective this way will make the change of magnification to be easier during using.

Clean the objective regularly, for lens is susceptible to dust.

When operating, use 10xmagnification objective to search and focus specimen firstly, then replace with higher magnification objective if necessary.

When replacing the objective, slowly turn the nosepiece until you hear "clicked", that means the objective is in place.

3.2.4 Mounting the filters

1. Open the condenser carriage①;
2. Place the required filter②into the condenser,then close the condenser carriage。

The filter of the standard outfit is green and baby blue.

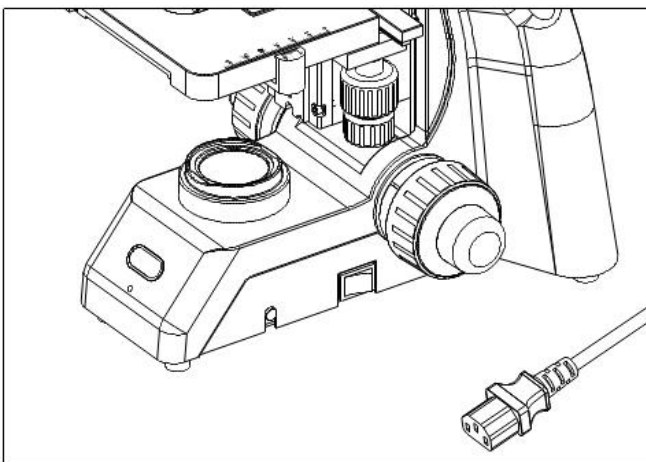


Figure 8

3.2.4 Connecting the Power Cord

The cable and cords are vulnerable when bent or twisted, never subject the power cord to excessive force.

1. Set the main switch①to "O" (off) state before connecting the power cord.。
 2. Insert the plugs②into the power jack③of the microscope safely.
 3. Plug the power cord④into the power supply receptacle⑤. Make sure the connection is well.
- Do use the supplied power cord all the time. If lost or damaged, select the same standard cord, please. A wide range of voltage ,like 100V~240V, is acceptable for this microscope.

3.2.5 Replacing the fuse Do remember to set the main switch①to the state "O" (OFF) and unplug the power cord before replacing the fuse. Rotate the fuse out of the holder③, replace with a new fuse, then rotate it back to the holder again.

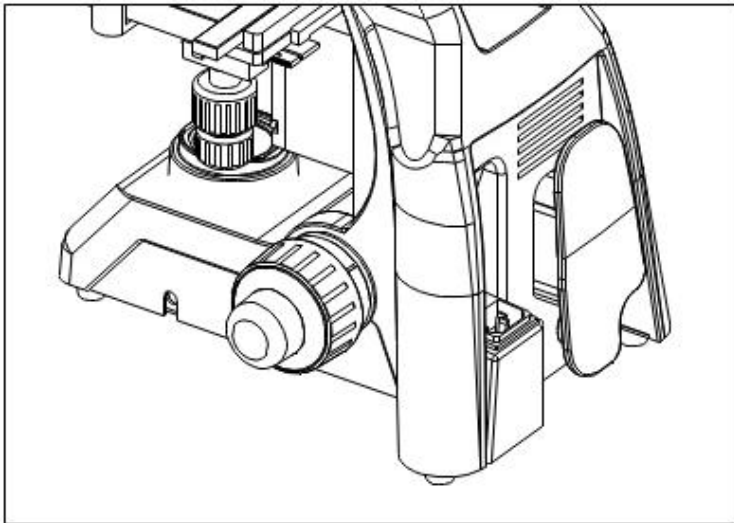


Figure 9

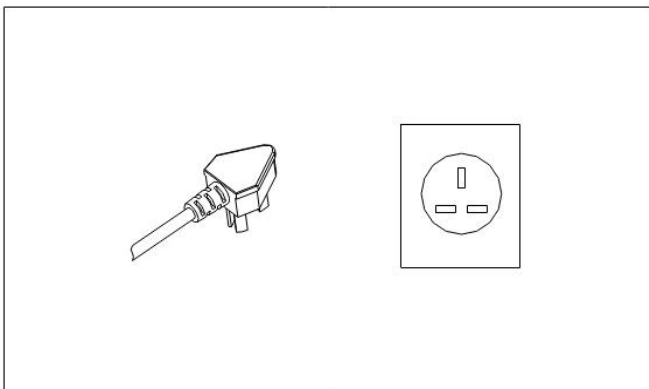


Figure 10

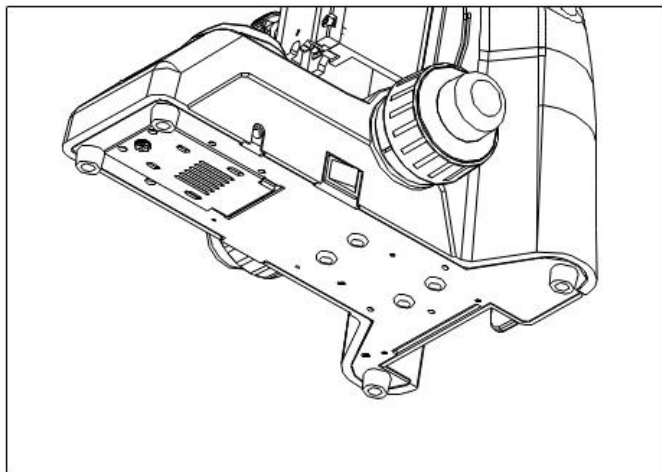


Figure 11

2.8 Installing and Replacing the bulb

Please use the specified halogen Lamp 6V20W.

1. Loosing the screw①.
2. Hold the bulb②after you wrap it with gauze or other protection materials, and then insert its pin deeply into the jack in the lamp holder.
3. During microscope using or soon after it is turned off, the bulb, the lamp house and nearby parts will be very hot and will cause serious burns. Please turn the main switch to "O" (off) and disconnect the power plug, and make sure the bulb, the lamp room and periphery are all cool. Then, you can do your replacing.

Insert the bulb gently. Squeezing too hard will damage the bulb.

To prevent reduced bulb life or cracking, do not touch the bulb with bare hands. If fingerprints are accidentally left on the bulb, wipe the bulb with a soft cloth.

4. Adjustment and operation

4.1 Adjustment set diagram

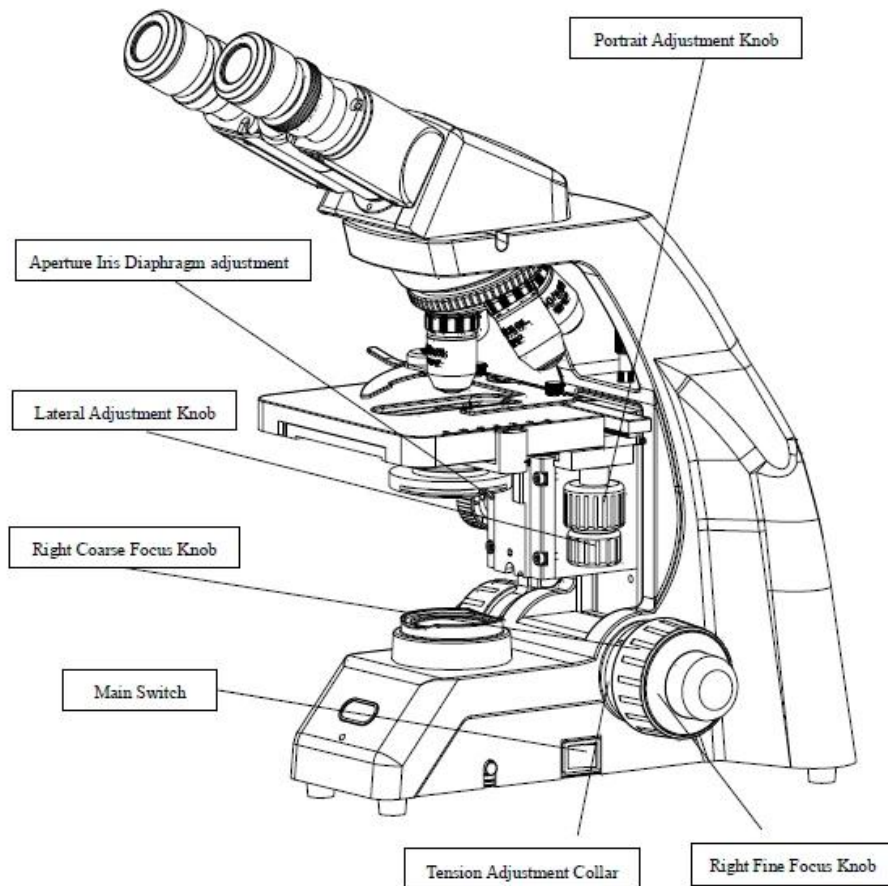


Figure 12

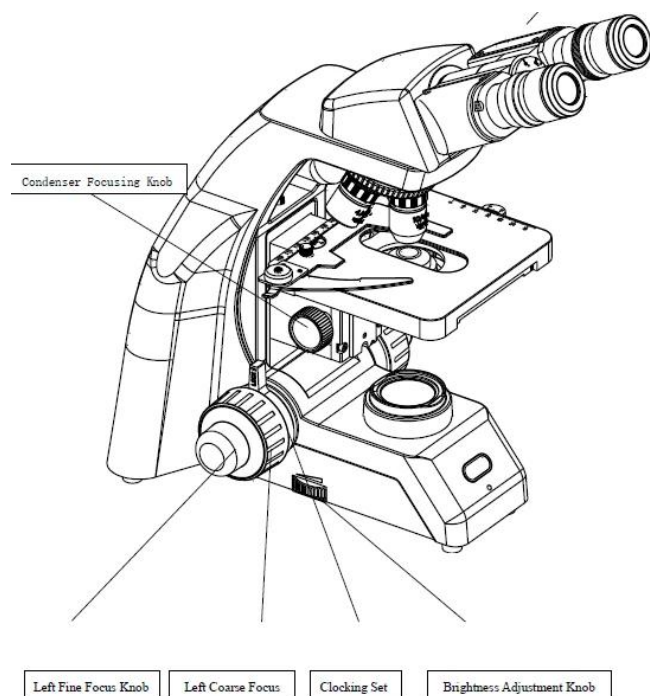


Figure 13

4.2 Adjustment and Operation

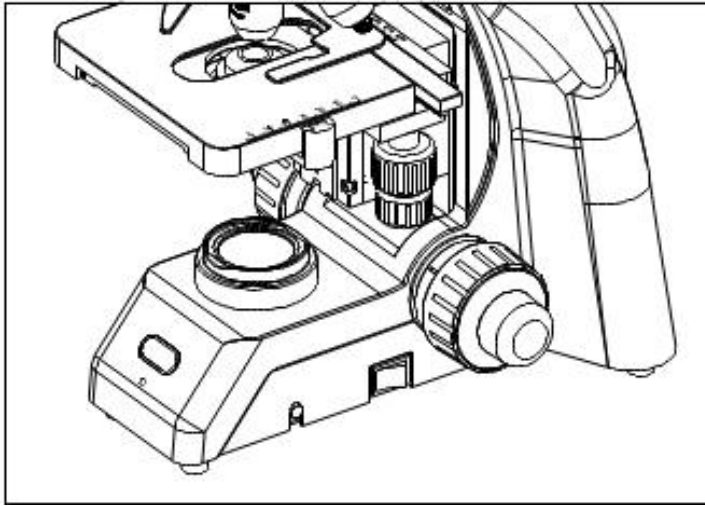


Figure 14

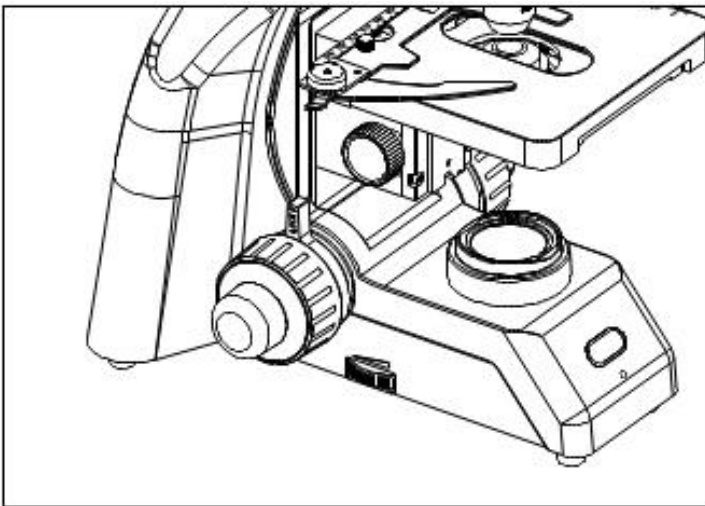


Figure 15

4.2.1 Brightness Adjustment

1. Connect the power cord and set the main switch④ to "-"state (ON).
2. Turning the brightness adjustment knob ②.

4.2.2 Placing the Specimen

1. Place the specimen the center of the mechanical stage and use the stage clips to clamp it.
2. Turn the portrait and lateral adjustment knob①of the mechanical ruler, move the specimen to the required position.

Be careful when changing the objective. If you finish the observation with the short working distance objective, and want to change another one, be careful of not letting the objective touch the specimen.

4.2.3 Focusing the Specimen

1. Focus the specimen with 10X objective. To avoid the objective touching the specimen during focusing,

you should raise the mechanical stage to let the specimen close to the objective at first, then slowly separate them to bring the specimen to focus. Turn the coarse focus knob① conversely to lower the specimen and search images in the 10xocular simultaneously, and then use the fine knob② to focus. After that, you can replace with other magnification objectives safely, and focus without the risk of damaging the specimen.

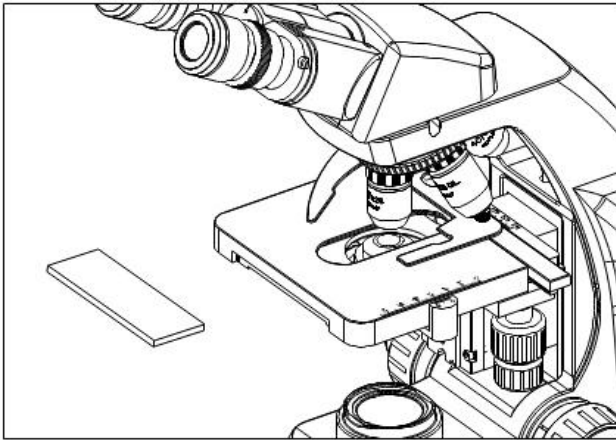


Figure 16

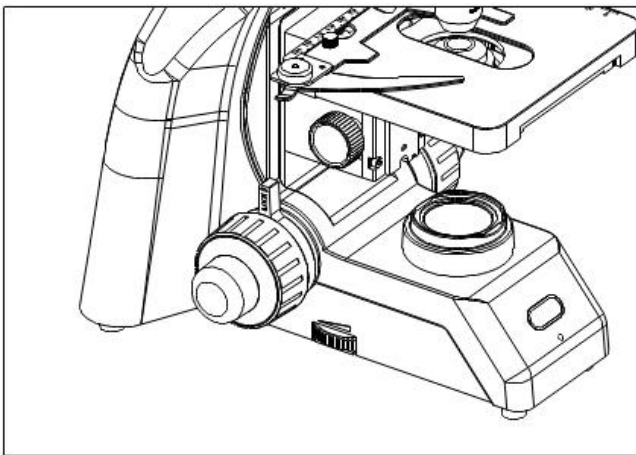


Figure 17

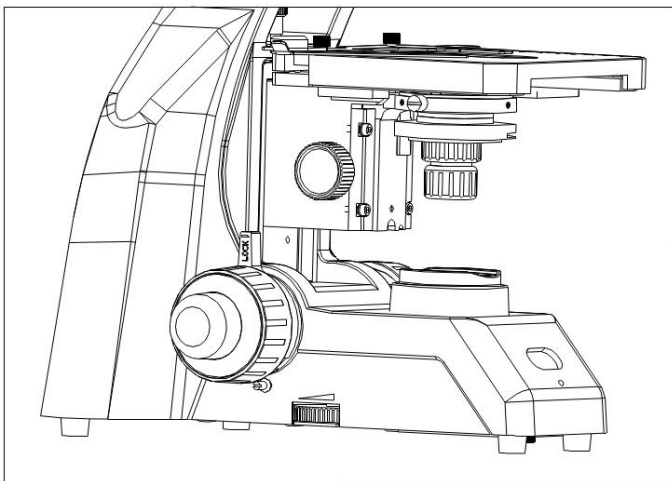


Figure 18

4.2.4 Condenser Adjustment

Turn the condenser focus knob①to move the condenser up and down. Raise the condenser when using the high magnification objective, and descend it when using the low magnification one.

1. Focus the specimen with 10x objective.
2. Adjust the condenser focus knob①to get a clear image of the field iris diaphragm.
3. Turn the condenser centering knobs②to center the image of field iris diaphragm in the field of view.

The condenser and the objective are coaxial. It has been adjusted before leaving factory, so the user needn't to adjust them by self

The highest position of the condenser has been adjusted too. It also needn't any user's operation. (The top surface of the condenser is 0.03mm-0.4mm lower than the stage top surface.)

4.2.5 Aperture

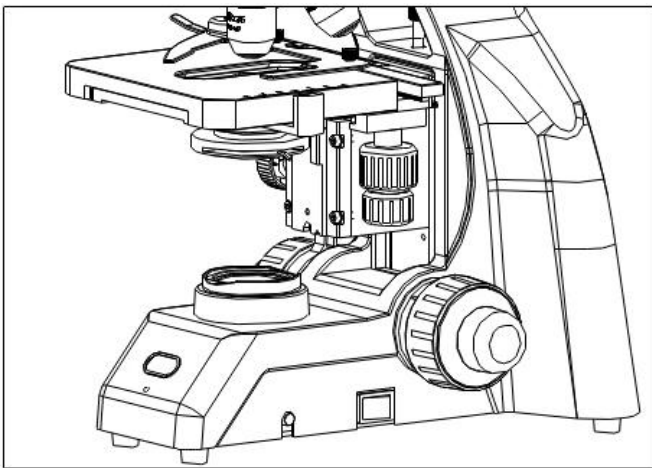


Figure 19

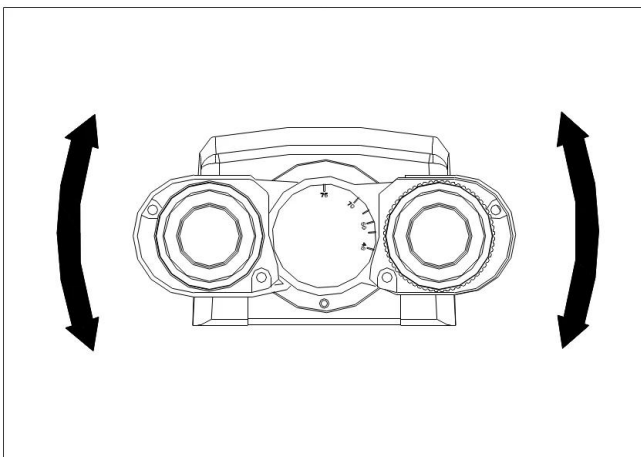


Figure 20

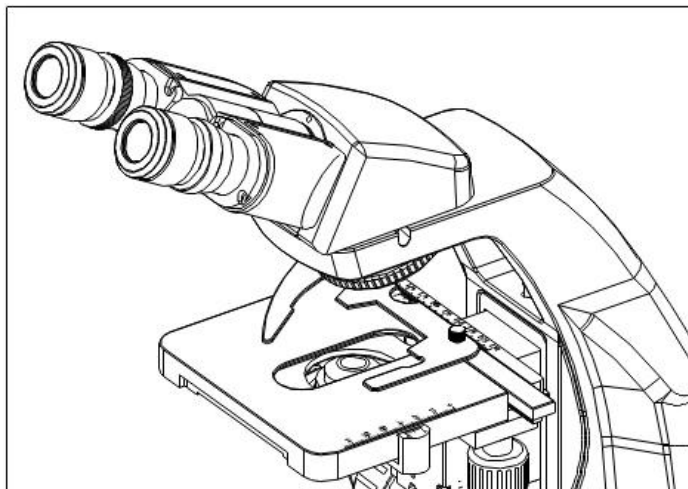


Figure 21

4.2.6 Adjusting the Interpupillary Distance

The interpupillary distance range: 55mm~75mm.

While looking through the eyepieces, move both eyepieces round until the left and right fields of view coincide completely.

4.2.7 Adjusting the Diopter

1. Turn the eyepiece① to adjust the diopter while looking through it.

The diopter range of the eyepiece is ± 5 diopter. The number aligned to the line on the viewing head is the diopter in use.

4.2.8 Adjusting the tension adjustment collar

Turn the tension adjustment collar① with your fingers. When the collar is turned in the direction of the arrow, the tension of the coarse adjustment knob② increases. Turning the collar in the opposite direction decreases the tension.

If the nosepiece descends on its own or if the specimen gets out of focus quickly even when it is brought into focus using the fine adjustment knob③, it means the tension of the coarse adjustment knob is too low. Turn the collar in the direction of the arrow to increase the tension.

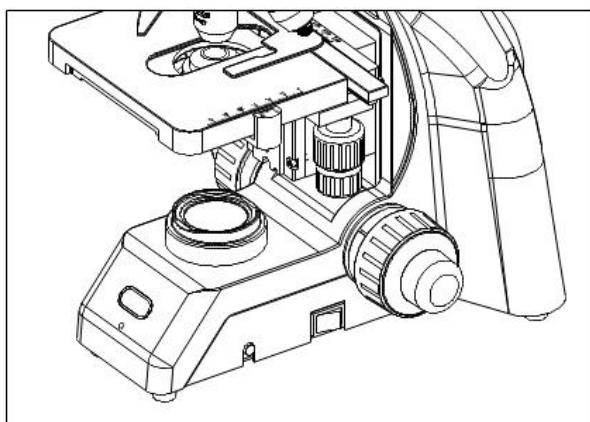


Figure 22

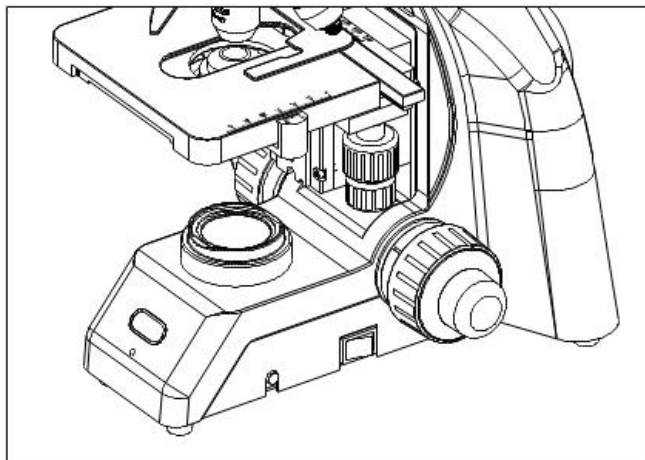


Figure 22

4.2.9 Adjusting the Interpupillary Distance

The microscope with ECO: The equipment can induct the front surroundings in 1m. At the moment, it light red and light source keeping working.

When people left and no object within the front 1m, the red light would flashed. After 15 minutes, light source stop working.

5. Technical Specifications

5.1 Main Specifications

Optical System	Limited optical system/Infinite optical system
Viewing Head	Seidentopf Binocular Head, 30°Inclined, Interpupillary Distance 48-75 mm
Eyepiece	WF 10X/18/WF 10X/20
Nosepiece	Backward Quadruple Nosepiece
Objective	Achromatic Objective/Infinite Semi-plan Objective 4x, 10x, 40x, 100x
Focusing	Coaxial Coarse and Fine Adjustment, Moving Range 0.002mm, Fine Division 20mm
Condenser	Abbe Condenser, NA1.25
Stage	Double Layers Mechanical Stage 160x140, Moving Range 74x50
	Timing Belt Stage 150x139mm, Moving Range 75x52mm
Illumination	1W S-LED /6V20W Halogen Lamp

Table 1

5.2 Eyepiece and objectives

5.2.1 Achromatic Objective

Magnification	Numerical Aperture(NA)	Focal Length(mm)	Focal distance(mm)	Working Length(mm)	Objective
4x	0.10	0.17	31.05	37.5	dry
10x	0.25	0.17	17.13	7.316	dry
40x	0.65	0.17	4.65	0.632	dry
100x	1.25	0.17	2.906	0.198	oil

Table 2

5.2.2 Infinite Semi-plan Objective

Magnification	Numerical Aperture(NA)	Focal Length(mm)	Focal distance(mm)	Working Length(mm)	Objective
4x	0.10	0.17	45	16.8	dry
10x	0.25	0.17	18	5.8	dry
40x	0.65	0.17	4.5	1.43	dry
100x	1.25	0.17	1.8	0.13	oil

Table 3

5.2.3 Eyepiece

Eyepiece	Magnification	Focal Length(mm)	Field of View(mm)
Wide field eyepiece	10x	24.95	Φ18
Wide field eyepiece	10x	25	Φ20

Table 4

5.2.4 Total Magnification

Eyepiece	10x	10x	10x	10x
Objective	4x	10x	40x	100x
Total Magnification	40x	100x	400x	1000x

Table 5

6. Outfit

Item	Specification	Quantity	Standard Outfit
			NE125
Main Body	Main Standard	1	•
	Main Standard(battery charging)	1	○
	Main Standard(ECO)	1	○
	Double Layers Mechanical Stage	1	•
	Condenser Holder	1	•
Viewing Head	Seidentopf Binocular Head	1	•
	Seidentopf Trinocular Head	1	○
Condenser	Abbe Condenser, NA=1.25	1	•
Nosepiece	Backward Quadruple Nosepiece	1	•
Illumination	1W SLED	1	•
	6V20W Halogen Lamp	1	○
	Blue, Green Filter	1 each	○
	Spare Fuse 50T250V 500mA	2	•
Eyepiece	10x Wide Field Plan Eyepiece	2	•
Objective	Achromatic Objective 4x	1	•
	Achromatic Objective 10x	1	•
	Achromatic Objective 40x	1	•
	Achromatic Objective 100x(Oil, Spring)	1	•

Table 6

Note: "•" Standard Outfit, "○" optional

Item	Specification	Quantity	Standard Outfit
			NE126
Main Body	Main Standard	1	•
	Main Standard(battery charging)	1	○
	Main Standard(ECO)	1	○
	Synchrone Belt Stage	1	•
	Condenser Holder	1	•

Viewing Head	Infinite Seidentopf Binocular Head	1	•
	Infinite Seidentopf Trinocular Head	1	○
Condenser	Abbe Condenser, NA=1.25	1	•
Nosepiece	Backward Quadruple Nosepiece	1	•
Illumination	1W SLED	1	•
	6V20W Halogen Lamp	1	○
	Blue, Green Filter	1 each	○
	Spare Fuse 50T250V 500mA	2	•
Eyepiece	10x Wide Field Plan Eyepiece	2	•
Objective	Infinite Semi-plan Objective 4x	1	•
	Infinite Semi-plan Objective 10x	1	•
	Infinite Semi-plan Objective 40x	1	•
	Infinite Semi-plan Objective 100x (Oil, Spring)	1	•

Table 7

Note: "•" Standard Outfit, "○" optional

7. Troubleshooting Guide

7.1 Optical System

TROUBLE	CAUSE	SOLUTION
1. The edge of the field of view is dark or the brightness is not uniform	The nosepiece is not in the located position (objective and light path not coaxial)	Locate the nosepiece properly where it clicks
	The surface of the lamp becomes black	Change a new lamp bulb
	A lens (the objective, condenser, eyepiece or collector) is dirty.	Clean it thoroughly
2. Dirt or dust is visible in the field of view	Dirt/dust on the specimen	Replace with a clean specimen
	Dirt/dust on the eyepieces	Clean them

3. Visibility is poor Image is not sharp; Contrast is poor; Details are indistinct	Specimen is not covered	Add cover glass on it
	The thickness of the cover glass is not suitable	Use standard cover glass with thickness of 0.17mm
	Specimen is placed reversely	Turn it over
	Dry objective has oil on it. (especially for 40X objectives)	Wipe the oil
	A lens (the objective, condenser, eyepiece or collector) is dirty.	Clean it
	Immersion oil is not used with the 100x objective	Use specified oil
	Air bubbles existed in the immersion oil	Eliminate the bubble
	The aperture iris diaphragm is stopped down too far	Adjust the aperture iris diaphragm properly
	Dirt or dust on the eyepiece	Clean it
4. One side of image is blurred	Condenser is not properly centered	Center the condenser with the centering screw
	The nosepiece is not properly engaged	Engage the nosepiece properly
	The specimen is not clamped	Clamp it with stage clips
5. The brightness is not enough	The aperture iris diaphragm is too small	Adjust it properly
	The condenser is too low	Adjust it properly
	A lens (the objective, condenser, eyepiece or collector) is dirty.	Clean it

Table 8

7.2 Mechanical System

TROUBLE	CAUSE	SOLUTION
1. Field of view of one eye does not match that of the other	Interpupillary distance is incorrect	Adjust interpupillary distance
2. Observation is tiring	The diopter is not proper	Adjust the diopter properly
	The brightness of the illumination is not proper for eyes	Adjust the lamp voltage

Table 9

7.3 Electrical System

TROUBLE	CAUSE	SOLUTION
1.The bulb cannot light	No power supply	Check the power cord connection
	The pin of the bulb doesn't insert properly	Insert the pin deeply
	The bulb broken	Replace with a new one
2. The bulb burns out suddenly	The bulb is not the specified one; The voltage is too high	Use the specified bulb; lower the voltage
3. the illumination is not bright enough	the voltage is too low	Raise the voltage
4. Image flicks	The bulb is about to burn out	Replace with a new one
	The bulb is not inserted deeply	Check its connection

Table 10



Biolab Scientific Ltd.

Trillium Executive Center, East Tower, 675 Cochrane Dr, Markham, Ontario L3R 0B8, Canada
Email: info@biolabscientific.com | Website: www.biolabscientific.com