



TUBE ROLLER / ROTATOR

BTUR-401

INDEX

1. FOREWORD	2
2. INSPECTION	2
3. SAFETY FEATURES	3
4. PACKING	4
5. STRUCTURE DESCRIPTION	4
6. OPERATING INSTRUCTIONS :	5
7. WARRANTY INFORMATION	6

1. FOREWORD

Our digital roller mixer , this is a simple and reliable device. We ask you to consider technical requirements of exploitation to insure pleasant and continuous usage of Machine.

2. INSPECTION

Please check the machine and accessories when you open the packing at first time . For all questions please contact manufacturer or local representative .

3. SAFETY FEATURES

Please read this manual carefully before operate this machine.

1. It must be use the original adapter.
2. The machine should be placed with a flat and steady surface for operating.
3. The machine which have 3D shaking action please cover the tube then prevent liquid samples overflowing .
5. Please do not moving when the machine is working.
6. The liquid samples loading can not over the heaviest loads.

4. PACKING

Digital roller mixer	1unit
Power cord	1unit
Operation manual	1unit

Table 1

5. STRUCTURE DESCRIPTION

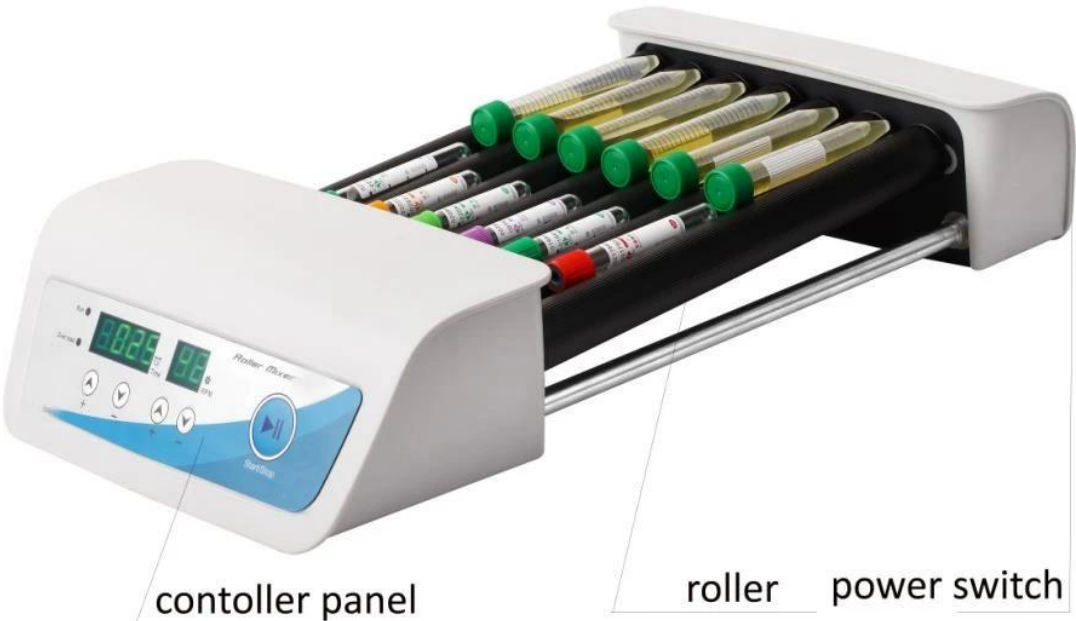


Figure 1

TECHNICAL SPECIFICATIONS:

VARIABLE SPEED REV/Min 1-99

SPEED SELECTION STEP REV/Min 1

Timing 0s-99 h

ENVIRONMENTAL TEMPERATURE, °C +2-50
AT THE TEMP, +20 °C): not more than 80% POWER ADAPTER INPUT : 110V-220V/50-60Hz POWER
ADAPTER OUTPUT: 12VDC,1.5A
POWER CONSUMPTION , no more than 20W

MODEL	DIMENSION (CM)	WEIGHT (KG)
BTUR-401	492*268*101	0.8

Table 2

6. OPERATING INSTRUCTIONS :



Figure 2

1. Press the power button after plug into a power . the display showing .
2. Press the "time adjustment button " setting the required operating time (Adjust the speed when machine operating is available)
Press the "Speed adjustment button " setting the required operating speed (Adjust the speed when machine operating is available)
3. Press the "START/STOP " button ,the machine start or stop operation .
4. If the overload light lighting or machine is stopping, please Press the "START/STOP" button and reduce the load or take out the tube which was stuck .

7. WARRANTY INFORMATION

Digital roller mixer is warranted for one year against defects in materials and workmanship

Should your digital roller mixer fail to function according to specifications at any time. Please contact the manufacturer or local representative immediately. The warranty will not cover defects caused by normal wear or by a breach of the manufacture's recommendations. Each Digital roller mixer is tested before shipping by the manufacturer. The quality assurance procedure guarantees that the roller mixer you have purchased is ready for use . Warranty repair could be performed only if the equipment is delivered in the original manufacturers packaging or equally safe packaging .Therefore please keep then transportation packaging after unpacking the device . If the above warranty requirements are disturbed ,repair charges are applied to the customer. For all further questions and maintenance please contact manufacturer or local representative .



Biolab Scientific Ltd.

Trillium Executive Center, East Tower, 675 Cochrane Dr, Markham, Ontario L3R 0B8, Canada
Email: info@biolabscientific.com | Website: www.biolabscientific.com