



# NITROGEN SAMPLE CONCENTRATOR

BCON-104

# INDEX

1. Introduction	4
.....	
1.1 Feathers	4
2. Specifications	5
.....	
2.1 Basic Parameters and Performance	5
3. Preparations	6
.....	
3.1 Structure Description	7
3.2 Instrument Installation	7
3.3 Installation steps	8
3.4 Installation of nitrogen flow rate adjusting valve	9
4. Operation panel	9
.....	
5. Operation Guide	10
.....	
5.1 Temperature and time set:	10
5.2 Operation and stop:	11
	12
5.3 Temperature error adjustment	12
5.4 Adjustment Methods:	13
6. Error Analysis and Recovery Processing	15
.....	
6.1 Error analysis and recovery processing:	16

## Foreword

Thank you for purchasing our Products: Nitrogen Sample Concentrator. This Manual for users contains function and operation of the Instrument. In order to use the instrument properly, please read this manual carefully before using the Instrument. Keep it for later use when you meet with difficulties.

### Opening Check

Please check the Instrument and Appendix with the packing list when you first open the instrument packing case. If you find there is something wrong with the Instrument and the Appendix, do contact the vendor or the producer.

## Safety Warnings and Guidelines

### Important Operation Information of The Security

Users should have an entire conception of how to use the instrument properly before operating it. Please read this operation manual carefully before using the instrument.

It is forbidden operating before read the operation manual. If you do not follow the instructions in the manual, the instrument may cause accidental injury during operation, and electric shock may occur. Read the guidelines and directions below and carry out the precautionary measure according to them.

### Security

To operation, maintenance and repair the instrument, please comply with the basic guidelines and the remarked warnings below. Otherwise, the instrument will suffer effect on the scheduled working life and also on the protection provided.



This product is a normal and an indoor using instrument, meet the standards of normal equipment of GB9706.1 Class I-B.



Before operation, read the manual carefully. These units are designed for using in the laboratory environments by who're knowledgeable in safe laboratory practices.



The operator should not open or repair the instrument by himself. Otherwise, the instrument will lose the qualification of repair guarantee or cause accidents. The company will repair the instrument based on warranty description.



A.C. power's grounding should be reliable to safeguard against an electric shock. The 3-pin plug supplied with thermo-shaker's power cable is a safety device that should be matched with a suitable grounded socket.



The temperature of metal block and top cover shell may be very high during the operation. There may be scald or boiling of the liquid. It is strictly prohibited any part of the body touching the instrument.



Close the test tube lid before put the tube into the block. Liquids may spill out in the block or


onto the device if the tube lid is opened, which will damage the block or the device.



Make sure the voltage is complying with the voltage required. Make sure the rated electrical outlet load no lower than the demand. Power line should be replaced with the same type if it any damage. Make sure there is nothing on the power line. Hold the jack when pull out the power line. Do not pull the power line. Do not put the power line in ambulatory place.



The instrument should be put in the place where of low temperature, little dust, no water, no sunshine or hard light, and of good aeration, no corrosively gas or strong disturbing magnetic field, and far away from central heating, camp stove and other hot resource. Do not put the instrument in wet and dusty place. The vent on the instrument is designed for aeration. Do not wall up or cover the vent. The distance between each device should be more than 100cm when there is more than one instrument.

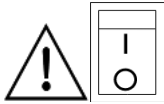

Main switch is on the rear of the device. Turn to “I” to power on the device, and Turn to “O” to power off the device.



Power off when operation finished. If long period do not use the instrument, pull off the connector plug, cover a cloth on the instrument to prevent from dust.



Pull the connector plug from the jack at once in the following case, and contact the vendor. There is some liquid flowing into the instrument; Drenched or fire burned; Abnormal operation such as abnormal sound or smell; Instrument dropping or outer shell damaged; The function has obviously changed

## Instruments Maintenance

The well in the block should be cleaned by the cloth stained with alcohol to assure good heat translation between the block and the test tube and no pollution. If there are smutches on the instrument, clean them with cloth.



- Power off when cleaning the instrument.  
Do not drop the clean fluid in the well when cleaning. Corrosive clean fluid is strongly prohibited.

## After-sales service

### 1- Warranty content

Within a month from the delivery date, the cause of failure due to material and manufacturing defects, the Company will be responsible for the guarantee replacement

Within 12 months from the delivery date, the cause of failure due to material and manufacturing defects. The company will prove to be defective instruments selected for repair or replacement

The product must be sent to the maintenance department of the company by the user. The freight will be borne by the user. The Company will bear the return shipping costs.

Over the warranty period, the Company will receive appropriate maintenance costs.

## 2) Warranty coverage

The above warranty is not suitable for improper maintenance, damage caused by unauthorized repair or modification.

# 1. Introduction

Nitrogen Sample Concentrator BCON-104 is mainly used for concentrating or preparing sample in batch. Such as drug screening, hormone analysis, liquid phase and mass spectrometry in the analysis of sample preparation. It works by blowing nitrogen, in the surface of sample which is being heated to accelerate evaporating and separating the solvent in the samples without oxygen.

## 1.1 Features

1. The heater causes the sample to be rapidly heated to the evaporation temperature, and at the same time, the gas is blown to the surface of the solution through the gas needle, which promotes the rapid evaporation of the solution and the concentration of the sample.
2. The height of the air chamber assembly is adjustable, and the standard air needle length is 150mm.
3. When the sample is concentrated, the liquid level position of the concentrated sample can be observed at any time.
4. Gas needle is controlled independently. Separately blow of each needle and flow regulating of each needle are available to avoid gas waste.
5. The entire equipment can be put into ventilation cabinet when the concentration sample in toxic solvents.
6. Built in overheat protection, automatic fault detection and fault beep alarm devices.
7. LED display immediate temperature and diminishing time. Operation is simple and convenient.

## 2. Specifications

1. The normal working condition-

The room temperature- 4°C ~ 45°C

The relative humidity- ≤70%

The using power- AC220V / AC110V ~ 50-60Hz

### 2.1 Basic Parameters and Performance

Model	BCON-104
Parameters	
Temperature Range	R.T.+5°C ~ 150°C
Temp. Setting Range	5°C ~ 150°C
Temperature Accuracy (40~100°C)	±0.5°C
Temperature Accuracy (100~150°C)	±1°C
Temperature Uniformity (100°C)	±0.5°C
Blocks Quantity	1 Standard Visible block
Time Range	1min~99h59min
Needle Plate Max. Lift Stroke	285mm
Gas-in Joint Outer Diameter	Φ7mm
Nitrogen Pressure	≤0.1MPa
Nitrogen Flow Rate	0~10L/min
Needle Length	150mm
Power	400W
Dimensions (mm)	260 x 220 x 525
Weight(kg)	5.8kgs

# 3.Preparations


This chapter introduces the instrument's mechanical structure, keyboard and each key's function and some preparations before power on. You should be familiar with this chapter before the instrument is first operated.

## 3.1 Structure Description



## 3.2 Instrument Installation

Assembly

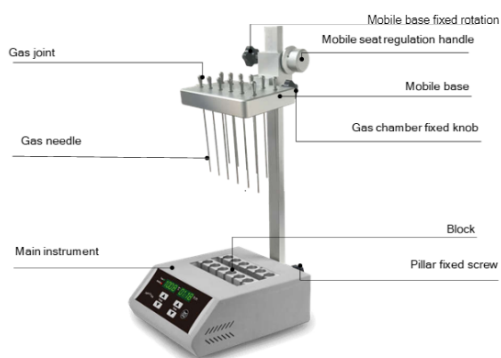


## 3.3 Installation steps

1. Put the dry bath on the bench smoothly and solidly
2. Fully insert the column into a square hole of the column holder behind the host, then screw knob into the column mounts, and finally tighten the knob to secure the uprights
3. Screw fixed knob of the mobile seat into the threaded hole on the left side of the mobile seat, then hold knob of mobile seat and counterclockwise rotate the hand wheel adjustment. When reach the desired position, we can tighten the knob to fix mobile seat. Clockwise rotating the hand wheel adjustment, the mobile seat will decline (as shown in the picture).

Note: When move the seat, two hands must be used in conjunction with operating mobile seat fixed knob and hand wheel adjustment

4. Screw fixed knob of the gas chamber into the threaded hole on the left side of the movable block, then insert column of gas chamber into the corresponding hole of the front movable block. Fully inserted column of gas chamber, then tighten fixed knob. (as shown in the picture)


### 3.4 Gas Control Valve (optional purchasing auxiliary)

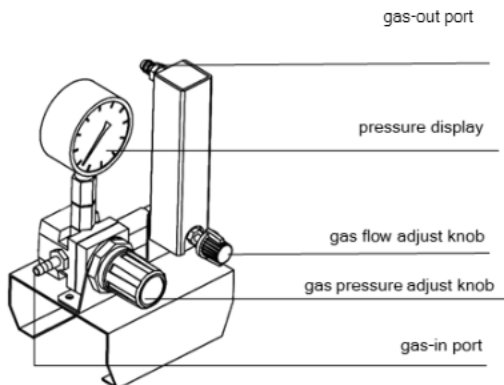

Pull out the gas pressure adjust knob, clockwise rotate the knob to increase pressure, anticlockwise rotate the knob to decrease pressure. Push back the knob to lock pressure. Clockwise rotate the gas flow adjust knob to increase flow rate, anticlockwise rotate the knob to decrease flow rate till shut off it.

## 3.4 Installation of nitrogen flow rate adjusting valve

The composition of nitrogen flow rate adjusting valve.

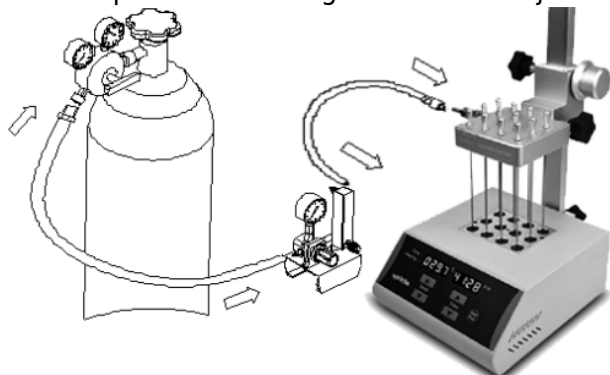


As the above picture, connect the outlet of the nitrogen flow regulator valve with outlet port of the host with a short hose (about 1.5m). Connect the inlet of nitrogen flow regulator valve with outlet port of the host with a long hose (about 3m)

Note- Slowly open the valve of the nitrogen bottle to control the outlet pressure between 0.1MPa to 0.2MPa, and then open the valve adjusting knob to keep pressure value at 0.02MPa. According to the number of holes, we can adjust the pressure appropriately (usually between 0.02MPa and 0.05MPa)



optional).

## 4. Operation panel

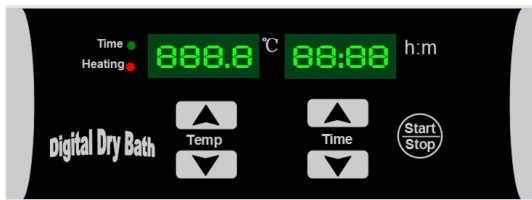

Button instruction:

- ▼ Decreasing button decrease the figure set
- ▲ Increasing button decrease the figure set

START/STOP Run/stop button - Run when you press the button once after setting the temperature and the time. Stop when you press this button for 2s.

## 5. Operation Guide

### 5.1 Temperature and time set:



The LED will display "8 " as the chart when the Instrument powers on and the Instrument goes into the initial state with the sound of "du...".




About 2s later, the figure 28.5 is the block's current temperature; 00:35 in the time display is the last set time.






Press ▲ or ▼ of the "Temp" key and keep off at once. Now, the value in the temperature display is the former setting temperature. As shown in the left drawing, the last digital of the setting temperature is flickering. If you want to set the temperature to 55.5°C, do as follows:



Keep pressing  when it reach 55.5°C, stop Pressing, it will auto save as 55.5°C  
Keep pressing the above keys for 2 seconds. We can modify the value fleetly. It is very convenient.



Press  or  of the “Time” key and keep off at once, the value in the time display is the former setting time. Shown in the left drawing is 00:35 (35minutes). At same time, the last digital is flickering.

If you want to change the time to 01:20, do as follows: Keep pressing , when it reach 01:20, stop Pressing, it will auto save as 01:20.

Note- If the time setting is 00:00, meaning the time of operation is forever, the Instrument run constantly in setting temperature.

## 5.2 Operation and stop:



Meanwhile, there is current temperature in the temperature display, and the radix point is flickering regularly during the rise process.  
When the temperature stops to rise, the flickering also stops. in the time display begins flickering, then the counting down starts.



When the timing is stopped, the operation is also stopped.

The buzzer alarms. Current temperature in the temperature display and **over** in time display mean “over”, the accomplishment of the operation.

After accomplishment of the operation, the instrument is in standby mode waiting for new declaration. Then pressing the keys near Temp or “Time” can reset the temperature or the time. Pressing START/STOP, it will operate according to last setting.

Press START/STOP for 2s during the operation, it will stop running. Press this button again, operation will continue.

## 5.3 Temperature error adjustment

The temperature of the Instrument has been adjusted before it is sold out. In case there is deviation between the actual temperature and the displayed temperature, you can do as follow for temperature calibration.

NOTICE:

1. The instrument has 2 calibration temperature points to ensure the veracity. It is linearly adjusted on 40°C and 100°C. The temperature accuracy will be within  $\pm 0.5^\circ\text{C}$  after temperature calibration. Both the environmental and the block temperature should be lower than 35° C when calibration.

## 5.4 Adjustment Methods:

- a- Power on the instrument, it enters into waiting interface. Make sure the current temperature in display is below 35°C. If the temperature is higher than 35°C, please wait until it down below 35°C.
- b- Inject olefin oil into one of block well, and then put a thermometer into this well (the precision of the thermometer should be 0.1°C and the temperature ball should be absolutely immersed into the olefin oil in the block well). Adiabatic material is needed on the block to separate it from the circumstance. (refer to below Figure A)

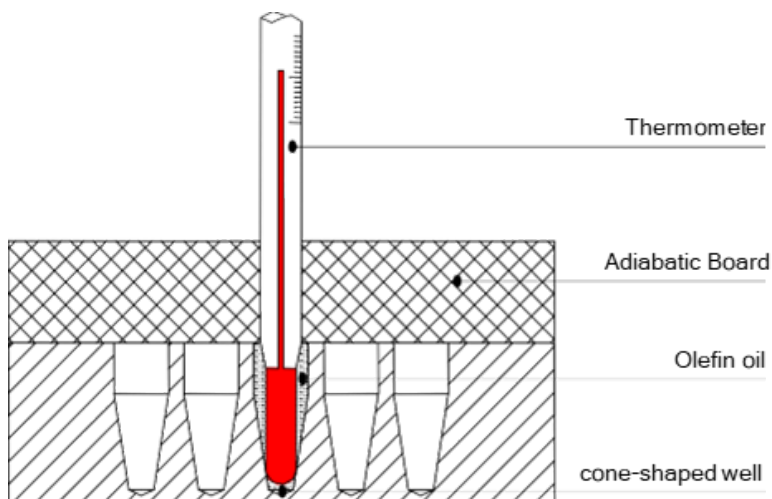




Figure A

NOTICE : To ensure the calibration precision, read the actual temperature value after the temperature reaches calibration points for at least 20 minutes.





Press  and  simultaneously to enter the temperature calibration interface as the right chart. LED display **Ad Jt** in the time display, the temperature display the current temperature and the program auto control the temperature to 40°C.



When the temperature reaches 40°C, the decimal digit begins to flicker. Wait for at least 20 minutes, after 20 minutes of constant temperature, read the measured temperature of the thermometer.

Notice- In order to ensure the temperature calibration accuracy, it is recommended that users read the temperature after 20 minutes of constant temperature  
Measured temperature.





If the thermometer reads 39.6°C, the actual temperature of thermometer is 39.6C  
press  and  of temperature to amend the temperature value to 39.6, Press "Start/Stop" to confirm.



Then the instrument automatically heats up to 100°C, and waits for the temperature calibration value to be input after a constant temperature of 100°C. It is also required to read the measured temperature of the thermometer after 20 minutes of constant temperature.

Notice! In order to ensure the temperature calibration accuracy, it is recommended that users read the temperature after 20 minutes of constant temperature  
Measured temperature





If the actual temperature of thermometer is 101.5°C, press  or  or of temperature to amend the temperature value to 101.5, Press "Start/Stop" to confirm.



After temperature have been adjusted, press  and  simultaneously to exit the temperature calibration and return to waiting interface.


Notice:

After temperature calibration, the temperature display is the same as the practical temperature of block.


Press  and  simultaneously during the calibration to exit the adjustment procedure, the calibration is invalid.

### The exchange of the block


a) Pull out the two screws which fix the block to the heating board with the screwdriver

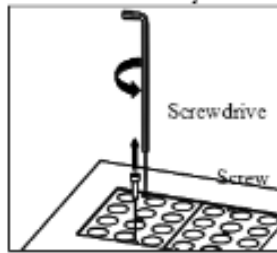



b) Fix the raiser in the center well of the block.

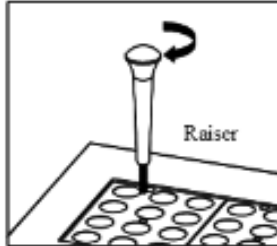

c) Pull out the raiser with the block.

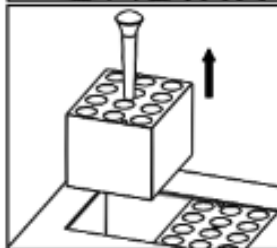

d) Screw the raiser out from the block, then fix the raiser to another block needed. Put the block in the instrument and fix it with screws.

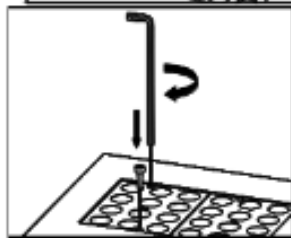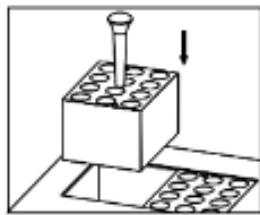

Notice: It is a little complicated to exchange the block, please be a bit patient

## 6. Error Analysis and Recovery Processing

### 6.1 Error analysis and recovery processing:


NO.	Fault phenomenon	Cause analysis	Recovery processing
1	No signals on the display when the instruction is powered on.	No power	Check the connection of power
		Broken fuse	Exchange fuse(250V 4A Φ5x20)
		Broken switch	Exchange the switch
		Others	Contact to the seller
2	The actual and displayed temperatures are quite different.	Broken sensor or loose contact of the block	Contact to the seller

3	" ERR " in the display with the alarm of "du..."	Broken sensor or the environmental temperature is below zero.	Contact to the seller
4	no heating of the block	Broken sensor.	Contact to the seller
		Solid state relay damage	
		Broken heater	
5	Key doesn't work	Key broken	Contact to the seller

Annex 1-

Wiring Diagram

(Below diagram is just for reference. It is subject to change without prior notice.)

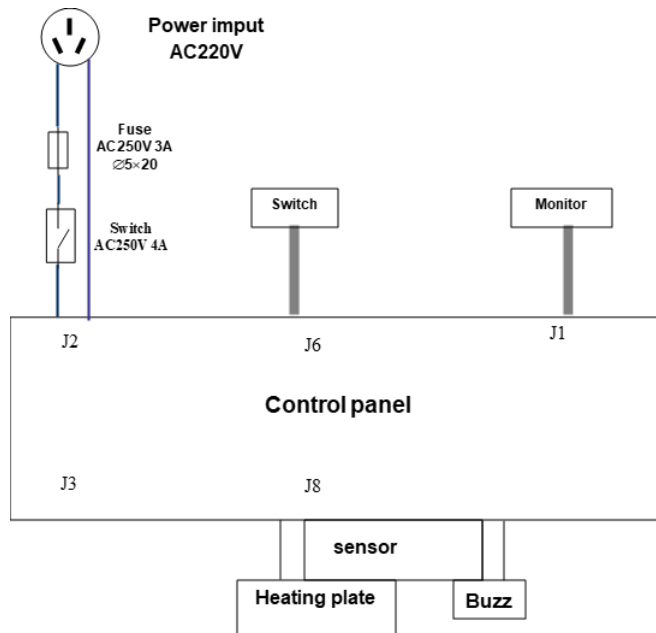

Biolab Scientific Ltd.

3660 Midland Avenue, Suite 300, Toronto, Ontario M1V 0B8, Canada  
 Email: info@biolabscientific.com | Website: www.biolabscientific.com